## CountAX TPC/CF

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Germ indicator for the determination of the total germ count (TPC) and for the coliformes screen (CF) from liquides and for surfaces.

---

### Gebrauchsanweisung Notice d'utilisation Instruction for use

**REF** AL0402

**Cont.** 10 Dip-Slides/Lames immergées/Dip slides  
10 Etiketten/Etiquettes/Labels

**Axon Lab AG**, CH-5405 Baden  
Tel. + 41 (56) 484 80 80  
Fax + 41 (56) 484 80 99  
info@axonlab.ch •  
www.axonlab.ch

# CountAX TPC/CF

### Intended Use


The CountAX Germ Indicator is used for the determination of the total germ count (TPC) and for coliform screen (CF) from liquids and on surfaces. The germ indicator can also be used for transporting samples.

### Package contents

- 10 Dip-slides
- 10 Labels

### Basic principle

The CountAX TPC/CF Germ Indicator includes 2 culture media: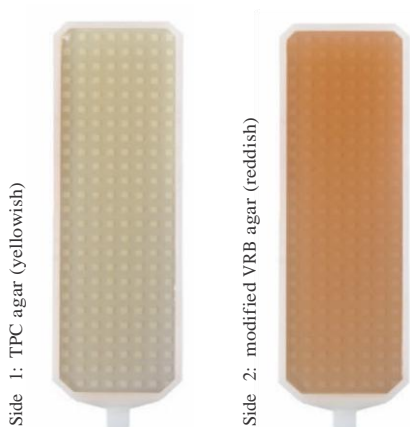

The TPC agar can be used to rapidly grow the most common bacteria and fungi.

The modified VRB agar side (crystal violet, neutral red bile agar with additional lactose) is used for the detection and enumeration of coliform Enterobacteriaceae. Lactose-positive Enterobacteriaceae are red/pink; lactose-negative Enterobacteriaceae are colourless.

### Composition

TPC agar	Modified VRB agar (CF)
Casein peptone	Peptone
Soy peptone	Yeast extract
Sodium chloride	Sodium chloride
Lecithin	Lactose
Polysorbate 80	Bile salts
Sodium thiosulfate	Crystal violet
L-Histidine	Neutral red
Agar	Agar

The culture media also contain substances that minimize the effect of disinfectants/detergents (surface tension).

### Storage

The CountAX TPC/CF will be stored at 15– 25°C. Temperature fluctuations and draughts must be avoided.

Do not freeze. Do not use the product after the expiration date (EXP) specified on the package.

Do not use the product if it is discolored or dried out, if the agar layers are detached from the plastic medium, or signs of bacterial/fungal growth. Do not touch the growing colonies as they might contain pathogenic germs.


### Sampling

Any contamination of the agar must be avoided. The agar surfaces must be fully placed on the analysis material. Once the sample has been used, screw the tube tightly.


## Performing the test

1. Unscrew the CountAX tube and remove the culture medium without touching the agar surfaces.

2. Contact inoculation

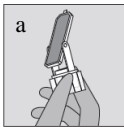
The flexible holder with the bracket enables easy handling (a). Solid surfaces can be tested by firmly pressing each side of the germ indicator against the relevant surface and holding it steadily for three to four seconds (b).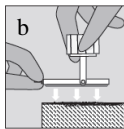

- 2a. Dipping process


Liquid samples are tested by dipping the germ indicator into the liquid for three or four seconds until the surfaces are completely covered (c). Use filter paper to wipe off any excess drops by the sample.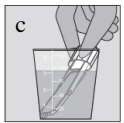

- 2b. Smearing process

Semi-solid materials or objects that are hardly accessible can be tested by smearing a sterile swab carefully for example over the frame of a limited area. If the object is dry, the swab will first be moistened with sterile water. A moistened swab can also be used to obtain samples from powders (e.g. spices) or viscous fluids.

After smearing the sample area, roll the swab carefully across the germ indicator's agar surfaces from left to right and bottom to top (d).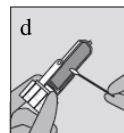

- 3.

3. Insert the culture medium in the tube and screw the tube tightly.
4. Complete the sample identification label and affix it to the tube.

5. Incubate the germ indicator tightly sealed for 1 day at 35 – 37°C or 2 days at 27 – 30°C or up to 5 days at 22°C.

## Transporting the inoculated CountAX tubes

The inoculated tubes can be transported at 15 – 25°C.


## Interpretation of the result

After incubation, remove the germ indicator from the tube, calculate the germ count (number of colony-forming units, CFU) and examine the color reactions. Lactose-positive Enterobacteriaceae are red/pink; Lactose-negative Enterobacteriaceae are colourless. The growth of gram-positive organisms is inhibited on the modified VRB agar.

The following limit values may be considered as a basis for evaluating the overall contamination level in food:

Good	< 10 <sup>5</sup> CFU/g
Permissible	10 <sup>5</sup> – 10 <sup>7</sup> CFU/g
Poor	> 10 <sup>7</sup> CFU/g

The limit values for cooked foods are lower than for raw foods. The presence of Enterobacteriaceae in cooked foods always indicates wrong handling or lack of hygiene.

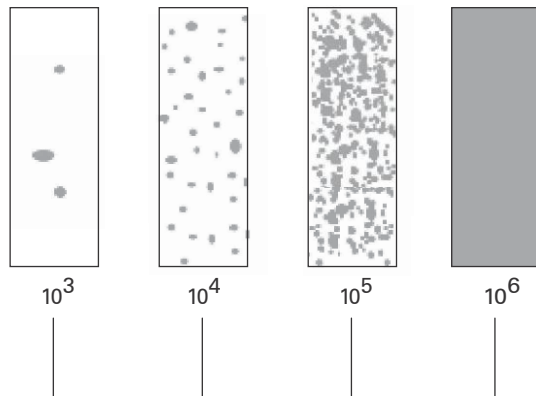It is difficult to specify CFU limit on surfaces since the criteria vary according to usage. However, a surface may be regarded as insufficiently cleaned if it shows 5 colonies/cm<sup>2</sup> after cleaning. This figure corresponds to about 45 colonies per slide on the TPC agar.

The following limit values may be considered as basis for evaluating the level of contamination:

Contact inoculation		
Clean	0	CFU/side
Contaminated	1–10	CFU/side
Highly contaminated	> 10	CFU/side

**Limitation of method**

When using the CountAX as a contact slide, the system is equally sensitive as the contact plate procedure.

The dipping and smearing processes have a detection limit of 1000 CFU/ml. The permissible TPC for drinking water is too low to be detected using CountAX slides.

Only results should be compared that were obtained by the same inoculation systems.

**Quality control**

Growth characteristics of the medium can be tested with the following strains:

	TPC	CF
Staphylococcus aureus ATCC	+	-
Enterococcus faecalis ATCC 29212	+	-
Escherichia coli ATCC 25922	+	+
Pseudomonas aeruginosa ATCC	+	+
Salmonella thyphimurium ATCC	+	+

(+) Growth (-) Partial or complete inhibition

**Waste disposal**

All laboratories are responsible for proper waste disposal in accordance with the legal requirements.

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